SVBONY

EN

DE

FR

IT

ES

RU

JP

SM403
STEREO MICROSCOPE
SERVICE MANUAL

Contents

A) Function	01
B) Features	01
C) Structure and Specification	01
D) Principle	03
E) How to Use	03
F) Maintenance and Care	23
G) Complete Sets	03
WARNING	04
GUARANTEE	29
EN(01-04)	
DE(05-08)	
FR(09-12)	
IT(13-16)	
ES(17-20)	
RU(21-24)	
JP(25-28)	

EU Importer: Germany Retevis Technology GmbH
Address: Uetzenacker 29,38176 wendeburg

A) Function

Use this stereo microscope for viewing, it has one pair of objective working simultaneously, so it is a binocular microscope with strong three-dimensional feeling. Its main function is as following:


1. As a research and dissection stool for zoology, botany, entomology, histology, mineralogy, archeology, geology, dermatology and so on.
2. Used for inspecting raw material and cotton fabric in textile industry.
3. As a welding assembly tool for mini-circuit boards, transistors, electronic components and so on in electronic industry, as well as their inspection tool.
4. For inspecting surface phenomena of various materials, such as crack formation, pore shape, corrosion and so on.
5. For inspecting surface quality of parts in the mechanical manufacturing. For inspecting the shape and size measurement of small precision parts, as well as their assembly tool.
6. For checking the surface quality of lens, prisms or other transparent substance, also for precision fine scale's quality checking.
7. For judging the true or false for the instruments and paper currency.
8. For jewelry and relics observation and identify.

B) Feature

1. This viewing head can rotate 360°, which greatly facilitates the user can observe at any location.
2. Long working distance makes greater operating space which is suitable for welding, assembly, dissection and so on.
3. This microscope is equipped with common auxiliary objective connector, can increase the magnification after coupled with auxiliary objective.
4. Rotate objective cover, select appropriate objective multiple.
5. When needed, can be screwed into auxiliary objective at the bottom of the lens cover.
6. If feel the torque too tight when using, rotate the focusing knob arduously; Or the torque too light, observing head auto-drop down, pls use wrench to adjust the torque of inside focusing knob to be appropriate.

C) Structure and Specification


1 . YJ-T101G Pole Type Stereo Microscope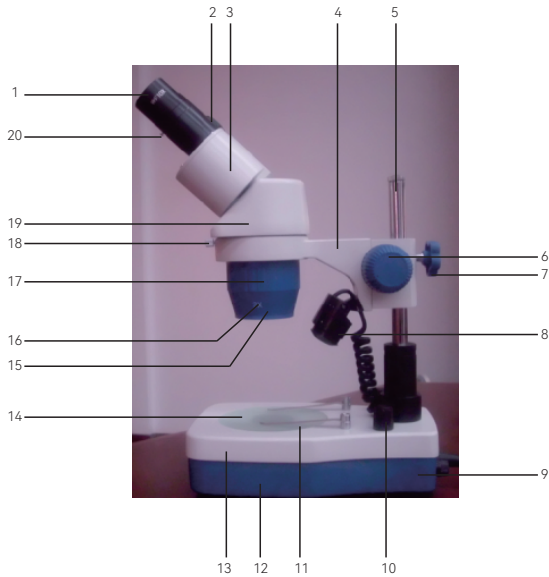

- | | |
|-------------------------------|-----------------------------|
| 1. Eyepiece | 11. Clips |
| 2. Diopter | 12. Lower Stage |
| 3. Prism Box | 13. Upper Stage |
| 4. Bracket | 14. Stage |
| 5. Pole | 15. Lens Dust Cover |
| 6. Focus Knob | 16. Objective Magnification |
| 7. Locking Screw | 17. Objective Cover |
| 8. LED Lamp | 18. Fixing Screw |
| 9. Switch | 19. Viewing Head |
| 10. Brightness Adjusting Knob | 20. Anti-theft Screw |

Category	Technical Specification
Eyepiece	Standard 1pair of WF10X (WF15X or WF20X optional)
Objective	Standard 2X, 4X (1X, 2X or 1X, 3X optional)
Viewing Head	Articulated binocular head, 45°inclined, 360°rotatable. Interpupillary distance 50-75mm adjustable, with a left diopter
Total Magnification	Objective Magnification X Eyepiece Magnification
Working Distance	100mm(The theoretical distance of lens from objects surface when the image is clear)
Base	Dimension 210X140X50mm, with clips
Stage	Diameter Ø 95mm, one frosted glass(using bottom light source), and one plastic black-white plate (using top light source)
Focusing	Rack and pinion focusing, with a focus torque handle wheel. Focusing Range: 40mm
Pole	Pole type, pole height 205mm. The moving range of head holder: 80mm
Illumination	Top and bottom high brightness LED lamp, brightness can be adjusted independently
Whole Size	210X140X360mm
N.W	2.5kgs
Package	Foam+ carton, package size 67X40X41cm(4sets/ctn)

2 . YJ-T101B Arm Type Stereo Microscope

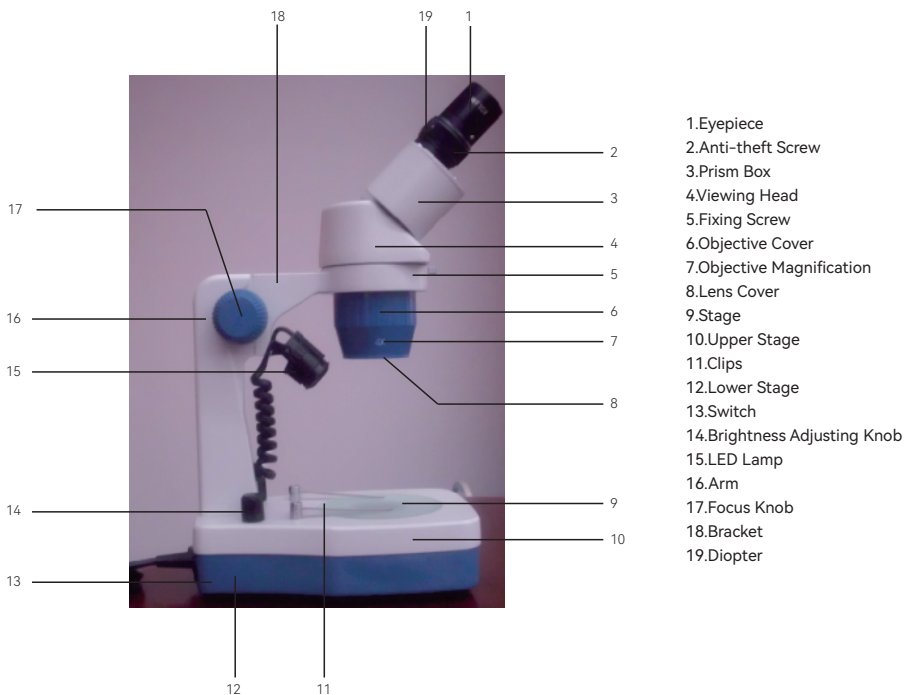Category	Technical Specification
Eyepiece	Standard 1pair of WF10X (WF15X or WF20X optional)
Objective	Standard 2X, 4X (1X, 2X or 1X, 3X optional)
Viewing Head	Articulated binocular head, 45°inclined, 360°rotatable。 Interpupillary distance 50-75mm adjustable, with a left diopter
Total Magnification	Objective Magnification X Eyepiece Magnification
Working Distance	100mm(The theoretical distance of lens from objects surface when the image is clear)
Base	Dimension 210X140X50mm, with clips
Stage	Diameter Ø 95mm, one frosted glass(using bottom light source), and one plastic black-white plate (using top light source)
Focusing	Rack and pinion focusing, with a focus torque handle wheel. Focusing Range: 40mm
Arm	Arm type, arm height 180mm.
Illumination	Top and bottom high brightness LED lamp, brightness can be adjusted independently
Whole Size	210X140X360mm
N.W	2.5kgs
Package	Foam+ carton, package size 67X40X41cm(4sets/ctn)

D) Principle

Using stereo microscope can get third dimension image, is due to can gather image from different directions by two objectives, then form different images on the eye's retina by two eyepieces. From the wide field eyepiece, can get a upright and third dimension image by binocular head. The left eyepiece tube with diopter ring, if the viewer's binocular degree are slightly different, you can adjust focusing knob so that right image to be clear, then rotate the left diopter so that the image to be clear. Binocular can be rotate in a certain angle to adjust the interpupillary distance.

This microscope has two pairs of different magnification objectives, can achieve different magnification's transform by turning the nosepiece (turn the objective cover), easy operation and fast, objective magnification are marked on the cover.

E) How to Use

1. Open the package and take out the microscope, insert the desired eyepiece, put on the rubber eyepiece cover and tighten the anti-theft screws(ensure the eyepiece free rotation)
2. Take off the objective cover, switch on. Select appropriate light source (top light or bottom light).
3. Put the objects in the centre of stag, then adjust the focusing knob. (When using YJ-T101G pole type stereo microscope, relax locking knob, then the observing head should be positioned in a appropriate height according to the objects' thickness, this distance should be about 70-80mm from the bottom of objective cover to the objects surface). Observe the right eyepiece to make the image clear(depending on the situation to select brightness). Adjust the left diopter again to make left and right images both clear.

F) Maintenance and Care

1. Whether the instrument in use or storage, should avoid dust, humidity, cold, heat or PH steam.
2. Chemicals shall not be placed near the instrument, even more in the microscope cabinet (except desiccant).
3. When not in use, the whole instrument should be covered by dust cover to keep it clean. If there is dust on the lens surface, don't wipe it by hand, only by brush.
4. If there is dust on the lens surface, can use clean old linen or cotton cloth dipped in a small amount of pure alcohol and ether mixture (1:1), or benzoline and xylene to wipe gently. Don't touch the lens surface by finger.
5. Due to time, grease on gear, rack, sliding groove surface and sliding cylinder will bond with dust into dirt or hardened, lead to rotating the coarse knob difficulty, can use xylene to clean, then add a little acid-free animal grease or acid-free vaseline, can't use other grease.
6. When microscope and eyepiece are not in use, should be covered with cloth or glass to prevent dust.
7. Do not put the plastic package a place your kids can reach. If not, the kids may put the plastic into the mouth and suffer from suffocating. Always supervise children using a optical device.
8. Store the product in a ventilated, dry, clean place or put it in a dry box.
9. Please do not disassemble or repair the product yourself. If you need to repair it, please contact our professional after-sales department.
10. Do not use non-waterproof products in wet environments, use waterproof products in wet environments, and dry them after use.

G) Complete Sets

- | | |
|--------------------------------|-------------------------|
| 1、 Microscope Body | 1set |
| 2、 Eyepiece WF10X | 1 pair (2pcs) |
| 3、 Eyepiece WF15X | 1 pair (2pcs, optional) |
| 4、 Eyepiece WF20X | 1 pair (2pcs, optional) |
| 5、 Rubber Eye Protective Cover | 2 pcs |
| 6、 Focus Torque Wrench | 1 pcs |

WARNING

Avoid Choking Hazard

WARNING:CHOKING HAZARD-Small Parts. Not suitable for children under 3 years old.

Authorized Accessories List

- Authorized Accessories meets the regulatory standards when used with the Svbonny accessories supplied or designated for the product.
- For a list of Svbonny-approved accessories for your item, visit the following website: <http://www.Svbonny.com>

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Post Code: _____ Email: _____

Remarks:

1. This guarantee card should be kept by the user, no replacement if lost.
2. Most new products carry a one-year manufacturer's warranty from the date of purchase.
3. The user can get warranty and after-sales service as below:
 - Contact the seller where you buy.
4. For warranty service, you will need to provide a receipt proof of purchase from the actual seller for verification

Exclusions from Warranty Coverage:

1. To any product damaged by accident.
2. In the event of misuse or abuse of the product or as a result of unauthorized alterations or repairs.
3. If the serial number has been altered, defaced, or removed.

Hong Kong Svbon Technology Co.,Ltd

Add: Unit B, 5th Floor, Gallo Commercial Building, 114-118
Lockhart Road, Wanchai, Hong Kong

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