
SVBONY

Wireless Microscope

User Manual

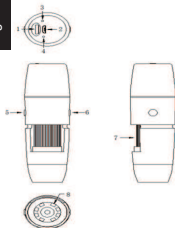SV602

1.The pictures are for reference only, please refer to the actual objects.

Part No.	Function
1	LED LED brightness level knob
2	Micro USB Micro USB charging interface
3	Indicator hole
4	Reset hole
5	On / off
6	Take a picture / video
7	Focal length adjusting roller
8	Light emitting diode

1.1 Functions and operation instructions of each component

LED brightness: under the working state, turn to adjust the brightness of LED light source.

Micro USB charging: when the power of the device is low, connect a charger to charge the device. It is not recommended to use the device during charging (it will reduce the service life of the device battery).

Indicator light: This hole indicates the working state of the equipment. Under the working state, the blue light is always on. When the USB is plugged in, the red light will be on, and when the USB is fully charged, the red light will be off.

Reset: when the equipment is stuck or cannot be shut down abnormally during operation, use a thin pointed object to stamp the reset button to shut down.

On / off: press this key for 2 seconds to turn on the machine, and press this key for 2 seconds to turn off the machine in the working state.

Photo / video recording: under the working state of the device, short press can take photos and automatically save photos, long press 2 seconds to enter the recording mode, then long press 2 seconds to exit the recording mode and save the recorded video, which can be viewed in the APP folder or in your iOS/Android device photo album.

Focus adjustment roller: under the working state of the equipment, rotate this roller to adjust the focus and focus the shooting object.

LED light source: brightness level knob can adjust its brightness.

2. Use Wifi digital microscope on your iOS/Android device

2.1 Download APP

IOS: Search for "HVViewing Image Analysis and Measurement Software" in App store App Store to download and install or scan the following two-dimensional code to select iOS version for installation.

Android: Scan the following two-dimensional code selection (Google) version (international users) or (China) version (Chinese users) to download and install or enter an address from the browser to download and install.

Download QR Code

Or enter the following address in the browser to download.

<https://console.cli.im/active/common/view/FVKkby>

2.2 Turn on the equipment

Press the device start key for 2 ~ 3 seconds, and the blue light will always be on, and the device will be turned on.

2.3 Connect Wifi digital microscope to your iOS/Android device

Open Wifi settings of iOS/Android devices, open Wifi, find a Wifi hotspot prefixed with HVCam-xxx and click Connect (no encryption), and return to the main interface of iOS/Android devices after successful connection.

2.4 Introduction and use of app interface

Open the APP and enter the APP main interface.

2.4.1 APP home page

Help: introduction of functions and operation pages.

Preview: Click to view the real-time screen of the equipment and operate the equipment.

File: Click to view the photos and video files already taken.

2.4.2 Preview interface

2.4.3 Key function description

Return: Click to return to the previous level.

Mirror left and right: Click to mirror the image of the device left and right.

Mirror up and down: Click to mirror up and down the device image.

Image flip: Click to cycle through the view.

Zoom out: Click to zoom out the screen (the default is the smallest for each opening).

Enlarge: Click to enlarge the picture (the picture is too small to use).

Center mark: click to mark the center of the picture with a cross.

Photography: Click to take photos and automatically save files.

Video recording: Click to record video / end video recording and automatically save the file.

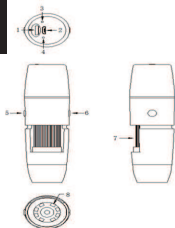
File: Click to view the photos or videos that have been taken.

2.4.4 Folder

Click on the folder to view photos or videos or delete photos or videos.

1. The pictures are for reference only, please refer to the actual objects.

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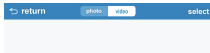
EU Importer:
Name: Germany Retevis Technology GmbH
Address: Uetzenacker 29,38176 wendeburg

EN:

 WARNING	<ul style="list-style-type: none">• Do not view the sun or other intense light with your products or even with the naked eye; Otherwise there may be permanent eye damage.• Kids need to be used under the guardianship of adults to avoid the unnecessary damage.• Do not put the plastic package a place your kids can reach. If not, the kids may put the plastic into the mouth and suffer from suffocating.
---	--

Care and Clean:

The item do not need routine maintenance other than making sure that the objective lenses and eyepieces are kept clean. Dirty objectives and/or eyepieces mean less light transmission and loss of brightness as well as unsharp images. Keep your optics clean! When not using your item, store those in the case provided.

Usage Notice:

1. Store the product in a ventilated, dry, clean place or put it in a dry box.
2. If there are dust and other dirt on the optical lens, use a professional cleaning kit to clean it. Do not wipe it directly with a cloth and a hard brush.
3. Please do not disassemble or repair the product yourself. If you need to repair it, please contact our professional after-sales department.
4. Do not use non-waterproof products in wet environments, use waterproof products in wet environments, and dry them after use.
5. When viewing indoors, please do not close the window and watch it through the window, which will reduce the product observation effect.

Avoid Choking Hazard

Small Parts. Not for children under 3 years.

Approved Accessories

- This item meets the regulatory standards when used with the Svbonny accessories supplied or designated for the product.
- For a list of Svbonny-approved accessories for your item, visit the following website: <http://www.Svbonny.com>

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Warranty Card	
Product Model	
Purchasing Date	
Defect Reason	
Dealer Name	
Telephone	
User's Name	
User's Address	
User's Email	

Remarks:

1. This guarantee card should be kept by the user, no replacement if lost.
2. Most new products carry a one-year manufacturer's warranty from the date of purchase.
3. The user can get warranty and after-sales service as below:
 - Contact the seller where you buy.
4. For warranty service, you will need to provide a receipt proof of purchase from the actual seller for verification

Exclusions from Warranty Coverage:

1. To any product damaged by accident.
2. In the event of misuse or abuse of the product or as a result of unauthorized alterations or repairs.
3. If the serial number has been altered, defaced, or removed.

MADE IN CHINA

Hong Kong Svbonny Technology Co.,Ltd

Add: Unit B, 5th Floor, Gallo Commercial Building, 114-118
Lockhart Road, Wanchai, Hong Kong

Facebook: [facebook.com/svbonny](https://www.facebook.com/svbonny)

E-mail: info@svbonny.com

Web: www.svbonny.com